



ROGYPAL AC-307-Z

Rinse cleaning compound

- **ROGYPAL AC 307-Z** is a specially developed rinse-cleaning compound for cleaning after using one of our alkaline cleaning concentrates, rinsing and neutralizing all types of glassware rinse in washing machines.
- Because **ROGYPAL AC-307-Z** is effective at rinsing, , no films remains behind which can be a breeding ground for bacteria.
- **Biodegradable.**
- **Phosphate-free.**

Where to use

Clear, liquid, odourless and colourless, acid rinse-cleaning compound for cleaning, rinsing and neutralizing of laboratory glassware and other equipment either manually or with a spray washer in laboratories, hospitals, institutions and industry.

Safe for steel, stainless steel, zinc, glass, porcelain and plastic.

Packing

Box of 2 x 5 kg
Can of 10 kg

How to use

Spray Washers:
0.2% **ROGYPAL AC-307Z** by volume. Temp .: 20-60 ° C.

overall:
0.5 to 1% by volume in water. Temp .: 20-80 ° C.

Properties

physical aspect	: clear liquid
pH at working concentration	: 2.5 ± 0.5
Odor	: mild
Solubility in water	: full

Safety

The safety information of this product is mentioned in our Safety Data Sheet (MSDS).
If using cleaning- and/or disinfectant products use always personal protective equipment

* Only to be used by professional users.